



## PRODUCT SPECIFICATION

### BETA SESSIONAL POLYCELLULOSE C-FOLDED POUCH WIPES STERILE



#### PRODUCT INFORMATION

##### PRODUCT DESCRIPTION

Beta Pre-saturated non-woven Polycellulose sterile wipes  
 Supplied C-folded in a pouch ready immediate use  
 Increased disinfectant and retention from re-sealable pouch  
 Specifically designed for use in Cleanrooms

##### PRODUCT FEATURES

Non-woven polycellulose wipe with sealed edges  
 Pre-saturated with Beta to ensure optimum disinfection  
 Supplied in a small pack size to minimise waste

##### MANUFACTURING

The product is pre-saturated in ISO 14644-1 Class 5 (in use) unidirectional airflow cabinets  
 The disinfectant is filtered through a 0.2 micron filter at the point of fill

##### PACKAGING

Supplied double bagged  
 All packaging is completed within an ISO 14644-1 Class 7 (in use) cleanroom

##### STERILISATION DETAILS

Sterilised by Gamma Irradiation (minimum 25kGy)

##### SHELF LIFE

Two years from date of manufacture

##### STORAGE CONDITIONS

Store in a cool dry place (< 40°C) away from direct sunlight and fluorescent light

## PRODUCT INFORMATION

### DIMENSIONS

Wipe Size: 23cm x 23cm (9" x 9")

### DISINFECTANT PERFORMANCE

| SUSPENSION TESTS                      |          |                            |                              |                    |            |                       |                        |
|---------------------------------------|----------|----------------------------|------------------------------|--------------------|------------|-----------------------|------------------------|
|                                       | Protocol |                            | Organism                     | Contact Time (min) | Conditions | Log reduction to pass | Log reduction achieved |
| Suspension Tests<br>Phase 2<br>Step 1 | EN 1276  | Bactericidal               | Salmonella typhimurium       | 1 min              | Dirty      | > 5.0                 | 6.30                   |
|                                       |          |                            | Escherichia coli             | 1 min              | Dirty      | > 5.0                 | 6.20                   |
|                                       |          |                            | Enterococcus faecalis        | 1 min              | Dirty      | > 5.0                 | 7.30                   |
|                                       |          |                            | Enterococcus hirae           | 1 min              | Dirty      | > 5.0                 | 6.60                   |
|                                       | EN 1650  | Fungicidal/<br>Yeasticidal | Aspergillus niger            | 2 mins             | Clean      | > 4.0                 | > 4.00                 |
|                                       |          |                            | Candida albicans             | 2 mins             | Clean      | > 4.0                 | > 4.00                 |
|                                       | EN 13704 | Sporicidal                 | Clostridium difficile spores | 2 mins             | Clean      | > 3.0                 | > 3.00                 |
| Suspension Tests<br>Phase 2<br>Step 2 | EN 13697 | Bactericidal               | Pseudomonas aeruginosa       | 5 mins             | Clean      | > 4.0                 | 5.70                   |
|                                       |          |                            | Escherichia coli             | 5 mins             | Clean      | > 4.0                 | > 6.56                 |
|                                       |          |                            | Staphylococcus aureus        | 5 mins             | Clean      | > 4.0                 | 6.02                   |
|                                       |          |                            | Enterococcus hirae           | 5 mins             | Clean      | > 4.0                 | 4.38                   |
|                                       |          | Fungicidal                 | Aspergillus niger            | 15 mins            | Dirty      | > 3.0                 | 4.08                   |

## ORDERING/PACKING INFORMATION

| Non-Sterile Code | Sterile Code    | Description   | Size        | Pack Qty           | Pack Qty | Case Qty |
|------------------|-----------------|---|-------------|--------------------|----------|----------|
|                  | WSBT02009 - A/I | Beta Sessional Polycellulose C-folded Pouch Wipes Sterile | 230 x 230mm | 15 wipes per Pouch | 15       | 10 x 15  |